



HCT-15 cells were transfected using the Neon™ transfection system and 0.5 µg of a plasmid encoding the EGFP. 24 hours post transfection, the cells were analyzed by light (A) and fluorescence microscopy (B).

#### Electroporation parameters

Pulse voltage (v)	Pulse width (ms)	Pulse number	Cell density (cells/ml)	Transfection efficiency	Viability	Tip type
1,130	40	2	5 x 10 <sup>6</sup>	88%	80%	10 µl
1,210	20	2	5 x 10 <sup>6</sup>	74%	80%	10 µl
1,150	40	2	4.6 x 10 <sup>7</sup>	75%	80%	10 µl

#### Cell information

Cell type/description	HCT-15 (gastrointestinal tract)/cell line
Characteristics/species	Adherent/human
Tissue origin	Colon
Media	Dulbecco's Modified Eagle's Medium (DMEM), with 4,500 mg/L D-glucose, L-glutamine, and 110 mg/L sodium pyruvate; 10% FBS
Morphology	Epithelial
Double time	—
Subculturing	Medium renewal: every 2 to 3 days
Culture condition	Temperature: 37°C; atmosphere: air, 95%; carbon dioxide (CO <sub>2</sub> ), 5%